

MTI GSL-1100 Tube Furnace

Operational manual ([pdf](#)) /
Temperature controller ([video](#))



Set single temperature:

- Turn on furnace and wait to see blinking "End"
-  to see "C 01" (initial temperature in °C)
Set using   and  to change digit
-  to see "t 01" (time in minutes to go to the next temperature)
Set to >30min to prevent overshooting
-  to see "C 02" - Set to desired temperature
-  to see "t 02" - Set to desired time (e.g., 600min)
-  to see "C 02" - Set to same temperature
-  to see "t 03" - Set to -121 to turn furnace off at end
-   simultaneously to show "stop"
-  to run program
- To pause at the current step, press  two seconds
- To view step and the time remaining 
- To stop the program, press  two seconds